milenia biotec

Milenia[®] Hybridetect
Detection of DNA and Protein

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Company Profile and Products

Milenia Biotec has been founded in the year 2000. The company is dedicated to the development, manufacturing, marketing and sales of rapid diagnostic tests.

The core product lines are tests for the diagnosis of allergies and inflammatory diseases.

Milenia Biotec GmbH is certified according to DIN EN ISO 9001:2000 and to the Quality Management Standard DIN EN ISO 13485:2000.

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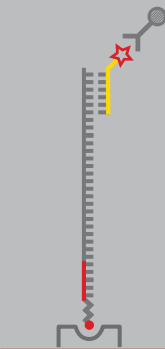
Principle of the Test (DNA-Detection) - A

1 DNA-Isolation

2 Addition of Primers + Nucleotides

3 PCR

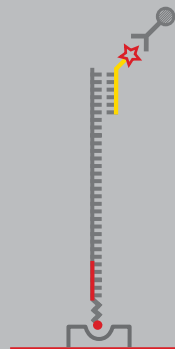
4 Addition of Probe

5 Hybridisation

6 Detection with Milenia Hybridetect

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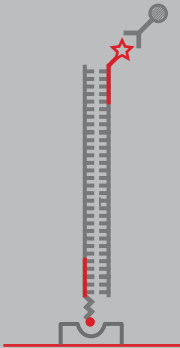
Principle of the Test (DNA-Detection) - B

1 DNA-Isolation

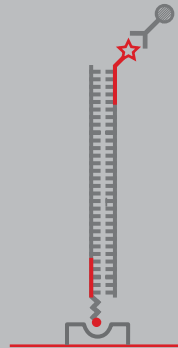
2 Addition of Primers + Nucleotides

3 PCR


4 Detection with Milenia Hybridetect

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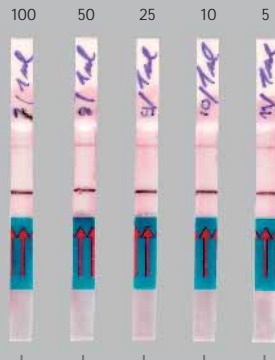
Characteristics of the Test and Example

Sensitivity: By using Milenia Hybridetect 5 pg of amplified DNA can be detected.

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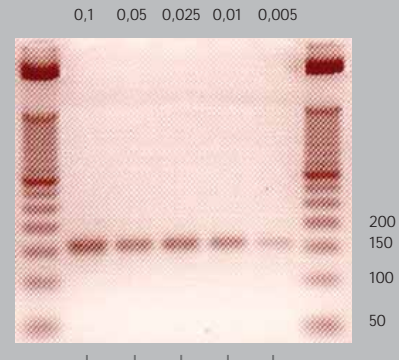
Compared to the previous Milenia Hybridetect test strip the sensitivity of the current strip (from lot 100) could be increased 10-fold.

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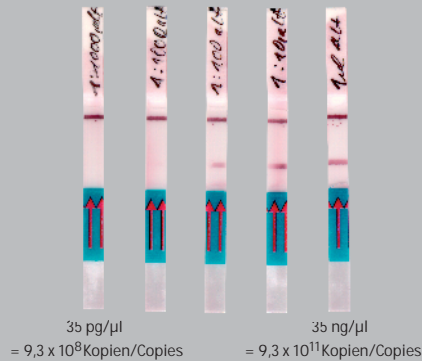Your Advantages:

- Time to Result: 15 minutes
- Robust Detection Method
- No Gel-Electrophoresis necessary
- Detection Limit: 5 pg DNA

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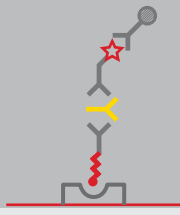
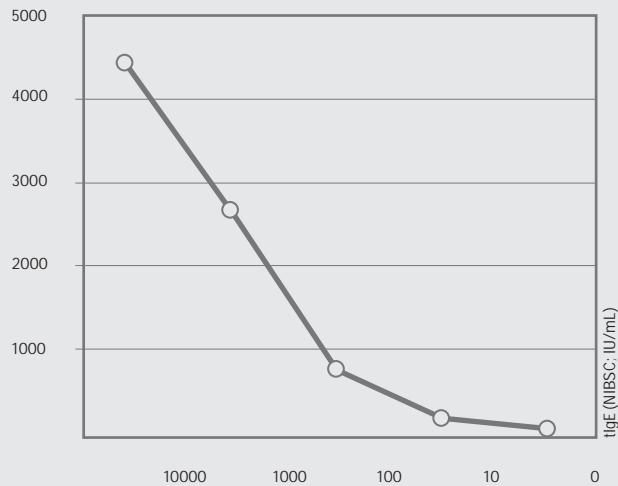
Characteristics of the Test and Example

Every immunoassay can be developed on the basis of the Milenia Hybridetect test strip, including the detection of proteins, small molecules and antibodies. The PicoScan instrument allows to report objective and (semi-) quantitative results. As an example an assay for the detection of total IgE is shown which is as sensitive as other immunoassays.

Order Information:

| Name | Order Code | Number Tests |
|-------------------------|------------|--------------|
| Milenia Hybridetect | MGHD 1 | 100 Tests |
| Milenia Hybridetect 2T* | MGHD 2 | 100 Tests |

* For the detection of 2 products on the same strip

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Allergy

Allergen specific IgE

- 29 Individual Allergens
- 2 Allergen Mixtures
 - Inhalation
 - Food

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
Total IgE

Immunology

- IL-6
- IL-6 Whole Blood
- IL-8
- LBP

Instrumentation

- PicoScan

milenia biotec

milenia biotec

Milenia Biotec GmbH

Hohe Straße 4-8 · D-61231 Bad Nauheim

Tel.: +49 6032 80 40-0 · Fax: +49 6032 80 40-80

E-mail: info@milenia-biotec.de

Internet: www.milenia-biotec.de

design: VISUV
fotos: Matthias Hoffmann Fotodesign

Allergie ▼ Autoimmun ▼ Immunologie ▼ Infektion ▼

HYB Milenia Biotec / A / 2007-07-30